

Public Lab


is an open community pursuing Do-It-Yourself environmental science.
This open hardware design was developed by Public Lab contributors; You are free to reproduce, share, & distribute with attribution, according to the CERN Open Hardware License 1.1.

[More instructions here](#)



Fold Here

make a diffraction grating from a DVD:
Cut a square from the outer edge of a DVD-R. The reflective part will peel off. We will use the clear part.

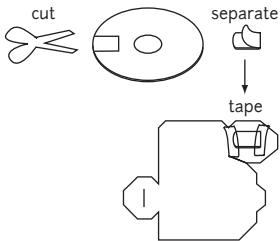

cut and fold:

Cut along the outer edge, then tape the DVD in place, oriented as shown.

Fold up or down as indicated by the dot and dashed lines. All labels should stay on the outside.

Tape flaps down onto the outside of the box.

Cut carefully and this card can be used to stencil more spectrometers.

